

IMPURITIES IN NITROGEN

Technique : Micro-GC

Column : Tamis

Carrier gas : Argon


Injector : Variable volume

Column temperature : NA

Injection time : NA

Column pressure : 35 psi

μTCD sensitivity : Standard

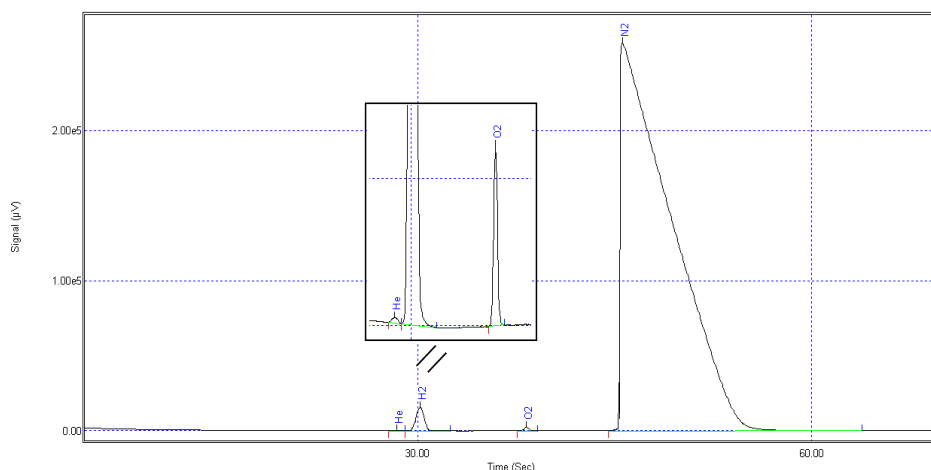

He	Hydrogen	Oxygen	Nitrogen
3.8 ppmV	792 ppmV	904 ppmV	99.83 %

The Micro-GC is a powerful solution that provides fast, accurate and easy-to-use solution to analyze your gas sample.

A wide range of capillary columns and MEMS injectors are available to optimize channel combinations inside the Micro-GC to obtain the desired separation. The μTCD is the high-performance universal detector used in such systems. A parallel analytical configuration of the channels or analytical modules will allow the sampling of a small amount of gas sample and simultaneous injection in all channels to develop up to 4 chromatograms and one cumulative report.

SRA Instruments is the ideal Partner with specific competences to develop with you turn key solutions based on optimized Micro-GC, accessories and dedicated software utility or report calculations for specific plant context.



More information : www.sra-instruments.com